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Contents

1. Identification data.....	9
2. General instructions.....	9
2.1. Symbols and means of representation.....	9
3. Safety.....	9
3.1. Grouped safety messages	9
3.2. Intended use	9
3.2.1. Liquids cabinet	9
3.2.2. Chemicals storage cabinet	9
3.3. Reasonably foreseeable misuse	9
3.3.1. Liquids cabinet	9
3.3.2. Chemicals storage cabinet	9
3.4. Personal protective equipment	9
3.5. Personnel qualifications.....	9
3.6. Duties of the operating company	10
4. Device overview	10
5. Transport	10
6. Setting up.....	10
6.1. Adjusting levelling base.....	10
6.2. Anchoring.....	11
6.2.1. Anchoring to the floor	11
6.2.2. Anchoring to the wall.....	11
7. Assembly.....	11
7.1. Chemicals storage cabinet.....	11
7.1.1. Containment tray.....	11
8. Maintenance	11
9. Cleaning	11
10. Spare parts.....	11
11. Technical data.....	11
12. Disposal.....	11

en

1. Identification data

Manufacturer	Hoffmann Supply Chain GmbH Franz-Hoffmann-Str. 3 90431 Nuremberg Germany
Brand	GARANT
Products	GridLine liquids cabinet GridLine chemicals storage cabinet, various versions
Version of the operating instructions	02 Translation of the original instruction manual
Date created	08/2020

2. General instructions

Read the instructions for use, follow them and keep them available for later reference.

2.1. SYMBOLS AND MEANS OF REPRESENTATION

Warning symbol	Meaning
 DANGER	Indicates a hazard which if not avoided will lead to death or serious injury.
 WARNING	Indicates a hazard which if not avoided may lead to death or serious injury.
 CAUTION	Indicates a hazard which if not avoided may lead to minor or moderate injury.
NOTICE	Indicates a hazard which if not avoided may lead to damage to property.
	Indicates useful tips and instructions together with information for efficient and problem-free operation.

3. Safety

3.1. GROUPED SAFETY MESSAGES

Chemical substances hazardous to the environment

Material and environmental damage in the event of improper storage of substances hazardous to the environment.

- » Do not exceed the maximum filling quantities.
- » Remove liquids that have leaked immediately and without leaving any residue.
- » Observe the material compatibility of the stored substances.
- » Only store chemicals in suitable containers and with proper labelling.
- » Keep doors closed.

Cabinet, roller cabinet or rack toppling

Risk of injury to hands, feet or body.

- » **It is essential the cabinet is secured against toppling.**
- » Wear safety shoes and safety gloves.
- » Do not store workpieces unsecured.
- » Do not open several drawers at once.
- » Keep cabinet shell cover clear of objects during transport.
- » Set down heavy objects slowly.
- » Ensure during movement and transport that all drawers and cupboard doors are closed and locked.
- » Comply with the maximum load capacity.

3.2. INTENDED USE

- For work in indoor industrial environments. For use on dry solid floors.
- Use only in a technically perfect and safe condition.
- Use exclusively approved Hoffmann Group accessories for modification or rebuilding.
- Keep doors' swing range unobstructed, keep doors and drawers closed.
- It is essential the cabinet is secured against toppling.
- Comply with the maximum load capacity.
- Secure stored containers against tipping over.

3.2.1. Liquids cabinet

- For storage of oil containers.
- When withdrawing liquids, position the oil tray under the container to be filled.
- Store a maximum of four containers of 25 litres in volume each or five containers of 10 litres in volume each per shelf.
- Oil container suitable for storing acid-free liquids, oils and greases with a flash point over 55°C.
- Only equip with pull-out shelf for liquids cabinets.

3.2.2. Chemicals storage cabinet

- For storing plant pesticides and liquids that are toxic and polluting to ground water of the GHS categories 1 to 4.
- For storing containers with a maximum filling quantity of 20 litres per shelf. Do not exceed total volume of 200 litres for all containers per shelf.
- Do not exceed the maximum load capacity of 100 kg per shelf.
- Only store substances that are compatible with the material of the containment tray.
- Only store compatible substances together.
- Maintain a minimum distance of 1 cm between the container and edge of the tray.
- Keep ventilation slots clear; do not cover.
- After maintaining the tray, have the manufacturer or specialised company carry out a leak test.
- Only equip with containment tray item no. 980096.

3.3. REASONABLY FORESEEABLE MISUSE

- Do not use in potentially explosive atmospheres.
- Do not make any unauthorised changes or modifications.
- Do not expose the work lamp to intense heat, direct sunlight or naked flames.
- Do not place objects on cabinet shell that will extend beyond its edge.
- Do not climb into or sit on cabinet, drawers or storage shelf.
- Do not place articles in the cabinet roughly.
- Do not store any foodstuffs.

3.3.1. Liquids cabinet

- Do not store any hazardous substances or chemicals.

3.3.2. Chemicals storage cabinet

- Do not store flammable liquids or aggressive substances.
- Do not use if containment tray has corrosion damage, cracks or leaks.
- Do not drill into containment tray or modify it in any way.
- No do expose it to water, precipitation or contamination.

3.4. PERSONAL PROTECTIVE EQUIPMENT

Comply with the national and regional regulations for safety and accident prevention. Select and provide protective work wear such as safety shoes and safety gloves appropriate for the risks associated with the intended activities. Wear suitable protective work wear when loading or withdrawing liquids. Whether the protective work wear is suitable or not depends of the hazardous substance used.

3.5. PERSONNEL QUALIFICATIONS

Specialists for mechanical work

Specialists in the sense of this documentation are persons who are familiar with assembly work, mechanical installation, commissioning, troubleshooting and maintenance of the products and who possess the following qualifications:

- Qualification / training in the field of mechanics as specified in the nationally applicable regulations.

Trained person

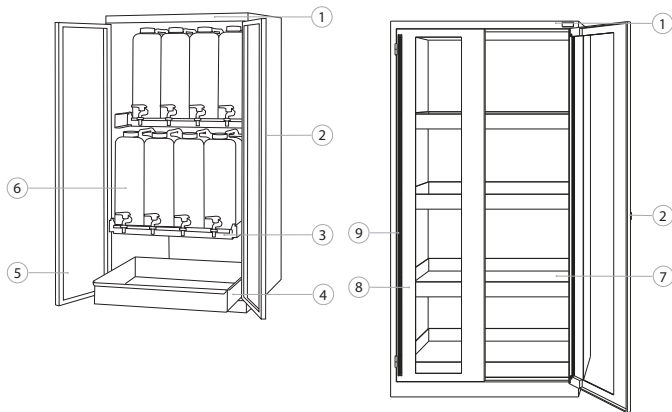
Trained persons in the sense of this documentation are persons who have been trained to perform work in the areas of transport, storage and operation.

3.6. DUTIES OF THE OPERATING COMPANY

The operating company must ensure that personnel who work on the product comply with the regulations and provisions together with the following instructions:

- National and regional regulations for safety, accident prevention and environmental protection regulations.
- No damaged products are assembled, installed or commissioned.
- The necessary protective equipment is provided.
- Observe the valid guidelines on storing hazardous substances.
- Observe directive concerning the requirements of containment trays made from steel (StawaR).
- Hazard assessment and preparation of a hazardous substance register by the operator are required.

4. Device overview

1	Cabinet shell	6	Oil container
2	Cylinder lock	7	Containment tray
3	Pull-out shelf	8	Viewing window swing door
4	Oil containment tray	9	Ventilation slot
5	Plain sheet metal swing door		

5. Transport

Immediately on receipt check the product for damage in transport. Damaged products must not be installed or commissioned. Pallet delivery. Transport on pallet using suitable lifting gear. Do not lift cabinet off pallet alone.

CAUTION

Risk of injury during transport

Risk of injury from improper transport to installation site.

- » Wear foot protection, safety gloves and safety helmet.
- » Do not transport alone.
- » Make sure the paths for movement and transport are clear.
- » If features, release parking brake on castors and apply before work.
- » Transport with drawers and doors closed, locked and additionally secured.

NOTICE

Improper transport

Damage to items of packaging.

- » Handle items of packaging in accordance with the symbols and instructions shown on them.
- » Do not drag them across the floor.
- » When items of packaging are being transported they must be upright, secured with ropes and protected against slipping.
- » Be aware of the centre of gravity.
- » Use the lifting points provided.
- » When putting them down, do so slowly and evenly.
- » Do not remove the transport packaging until immediately before erection.
- » Once the device has been erected at the place of erection, remove the transport restraints.

6. Setting up

- Set up on a solid, level floor suitable for the maximum load capacity of the cabinet.
- Use a spirit level to check horizontal levelling of the cabinet. If necessary adjust levelling base [► Page 10].
- All shelves and containment trays must be horizontal and level and protected against rain.

CAUTION

Chemical substances hazardous to the environment

Material and environmental damage in the event of improper storage of substances hazardous to the environment.

- » Do not exceed the maximum filling quantities.
- » Remove liquids that have leaked immediately and without leaving any residue.
- » Observe the material compatibility of the stored substances.
- » Only store chemicals in suitable containers and with proper labelling.
- » Keep doors closed.

NOTICE

Leaking liquids

Environmental damage and material damage.

- » When loading or unloading the cabinet, put the containers in intermediate storage properly.
- » Check containment tray and pull-out shelf for correct mounting and secure seating before loading.

6.1. ADJUSTING LEVELLING BASE

① Applies to: Chemicals storage cabinet.

NOTICE

Limited load-bearing capacity

Damage from overloading adjustable base.

- » When using the adjustable base, note load-bearing capacity reduced to 500 kg.
- » Before adjusting the adjustable base, completely clear out the cabinet.
- » Do not screw the adjustable base's thread all the way in. Keep 0.5 cm of thread protruding.

✓ Cabinet is completely empty.

1. Take lower storage shelf out of cabinet by lifting it at the holes on the sides.

CAUTION! Risk of cuts. Wear safety gloves.

2. Adjust four levelling feet inside the cabinet using a 5-mm hexagon L-wrench.
3. Use a spirit level to check horizontal levelling.
4. Repeat procedure until cabinet is perfectly level.

6.2. ANCHORING

① *Employ at least two persons when anchoring the cabinet.*

⚠ CAUTION

Tipping cabinet

Crush hazard for hands, feet and body.

- » **It is essential the cabinet is secured against toppling.**
- » Do not pull out multiple shelves at the same time.
- » Set down heavy objects slowly.
- » Comply with the maximum load capacity.

NOTICE

Wiring and pipes in wall

Damage to wiring and pipes.

- » Check wall for sufficient load-bearing capacity before drilling.
- » Use suitable drilling tools and materials.
- » Use suitable protective equipment.
- » Have trained personnel drill holes.
- » Do not drill holes close to wiring or pipes.

6.2.1. Anchoring to the floor

① *Applies to: Liquids and chemicals storage cabinet.*

📷 B

- ✓ Wall plugs and screws chosen to match the nature of the floor.
 - ✓ Floor has sufficient load-bearing capacity.
 - ✓ Adjusted levelling base. ▶ Page 10
1. Take lower storage shelf out of cabinet by lifting it at the holes on the sides.

CAUTION! Risk of cuts. Wear safety gloves.

2. Mark drilling spots through the holes in the cabinet floor and move the cabinet out of the way.
3. Drill the holes in the floor. Insert the plugs and bolt the cabinet to the floor.

6.2.2. Anchoring to the wall

① *Applies to: Chemicals storage cabinet.*

📷 C

- ✓ Wall plugs and screws chosen to match wall.
 - ✓ Floor has sufficient load-bearing capacity.
1. Mark drilling spots through the holes in the cabinet's rear and move cabinet out of the way.
 2. Drill holes in wall, insert wall plugs and screw cabinet to wall.
7. Assembly

⚠ CAUTION

Assembly work

Risk of crushing feet or hands when assembling and when closing drawers.

- » Wear safety shoes and safety gloves.
- » Ensure safe footing on clean, level surface.
- » Do not assemble by yourself.

7.1. CHEMICALS STORAGE CABINET

7.1.1. Containment tray

📷 D

① *Protective layer must not be damaged. In the event of damage, immediately remedy the fault or replace containment tray.*

1. Insert four shelf carriers per containment tray at an identical height in socket strip.
2. Place the containment tray on shelf carrier.

8. Maintenance

Interval	Action
Before the start of each shift	Check for externally evident damage, defects and for good condition. Check the edges and guides for signs of wear. <ul style="list-style-type: none"> ■ If defective, immediately block access and prevent use. ■ Have the damage repaired immediately.
Occasionally	Grease any locks and wheels with multi-function oil.
Weekly	Check for containers that have leaked. Remove liquids that have leaked from trays without leaving any residue.
Every two years	Check containment tray and underside of containment tray for visible damage, corrosion, leaks and damaged weld seams. Log result.

After maintaining the tray, have the manufacturer or specialised company carry out a leak test.

9. Cleaning

Do not use chemical, alcoholic, abrasive or solvent-based cleaning agents. Wipe down powder-coated surfaces (such as body, metal panelling, drawers, storage shelves) with lukewarm, dampened cloth. Clean polycarbonate windows with window cleaner and a soft, clean, non-abrasive cloth.

10. Spare parts

Order original spare parts from the Hoffmann Group Customer Service.

11. Technical data

	Load capacity	Drip catch volume
Pull-out shelf for liquids cabinet	200 kg	-
Oil containment tray for liquids cabinet	-	40 l
Containment tray for chemicals storage cabinet	100 kg	20 l
Overall load capacity of the chemicals storage cabinet	1000 kg	-

12. Disposal

Comply with the national and regional environmental protection and disposal regulations for correct disposal or recycling. Segregate items into metals, non-metals, composite materials and consumables and dispose of them responsibly.

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